



Sample Protein molecular weight standards Kit (Mw-GF-1000 KIT)

- | | |
|--|----------------|
| 1. Carbonic Anhydrase(Bovine Erythrocytes) | 1.5mg/mL 100μL |
| 2. Bovine serum albumin, BSA | 1.5mg/mL 100μL |
| 3. Alcohol Dehydrogenase, (Yeast) | 1.5mg/mL 100μL |
| 4. β-Amylase,(Sweet Potato) | 1.5mg/mL 100μL |
| 5. Thyroglobulin,(Bovine) | 1.5mg/mL 100μL |

Column	: Shodex PROTEIN KW-804 (8.0mmID*300mm) (Fig. 1)
	Shodex Asahipak GS-620 HQ (7.5mmID*300mm) (Fig. 2)
Eluent	: 0.1M Phosphate buffer(pH7.0) + 0.3M NaCl
Flow rate	: 1.0mL/min (Fig. 1)
	0.6mL/min (Fig. 2)
Detector	: MALS (Multi angle laser light scattering), Shodex RI
Column temp.	: 26°C